


Instructions Manual

Focusing Stand

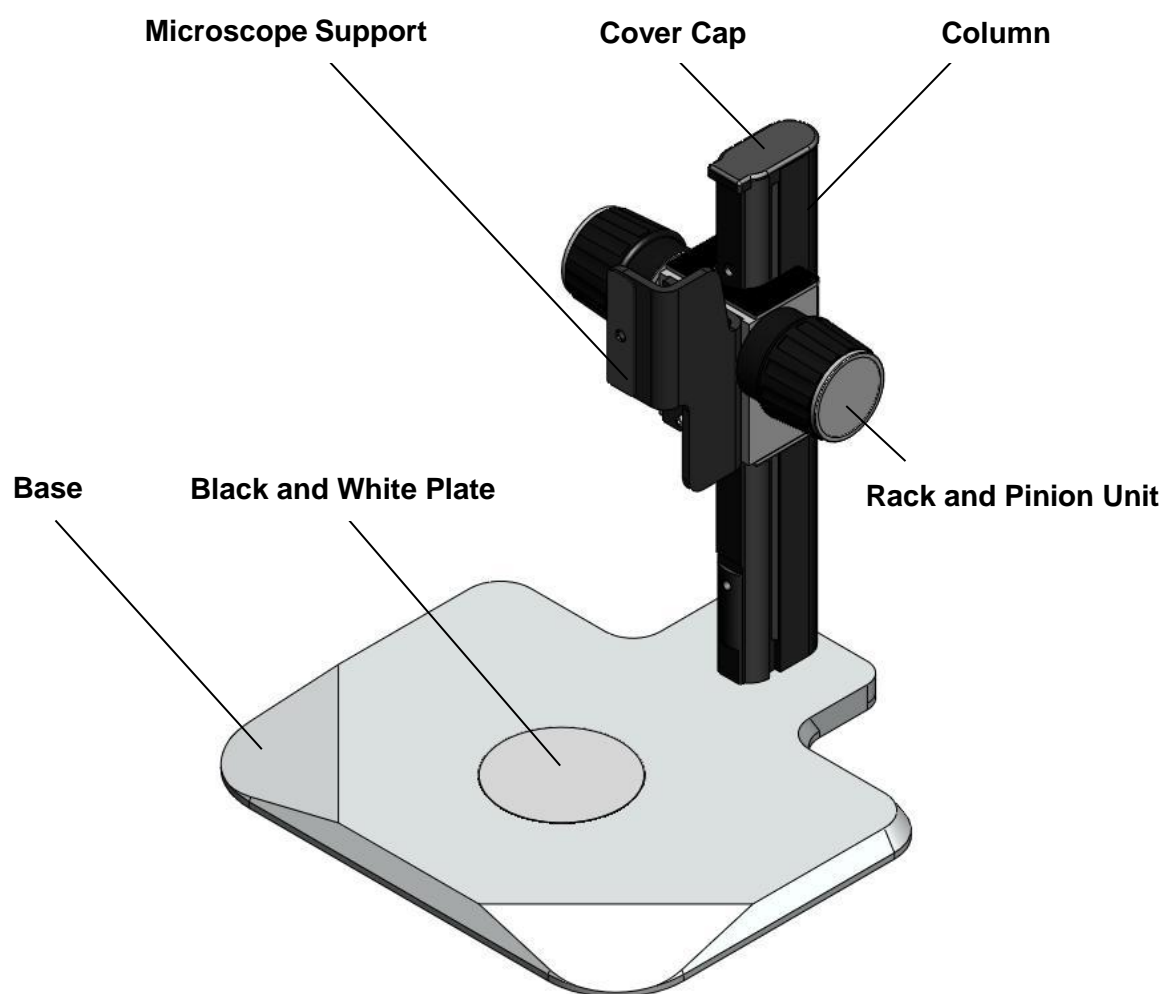


This manual describes how to install the Focusing Stand


Installation

Please follow the installation instructions below.


1. Place the Base on an even table by the side, use the included allen key to fasten the four screws on the bottom of the Column.

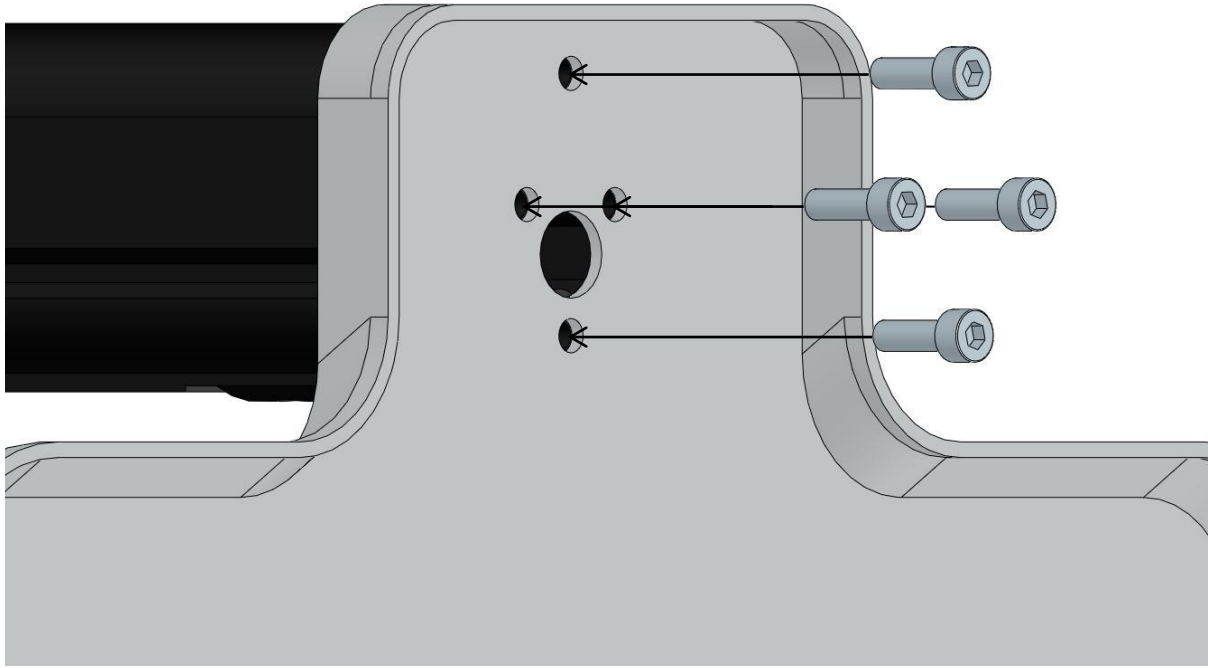

2. Install the Microscope Support Adapter on the Rack and Pinion Unit, and tighten the three screws.

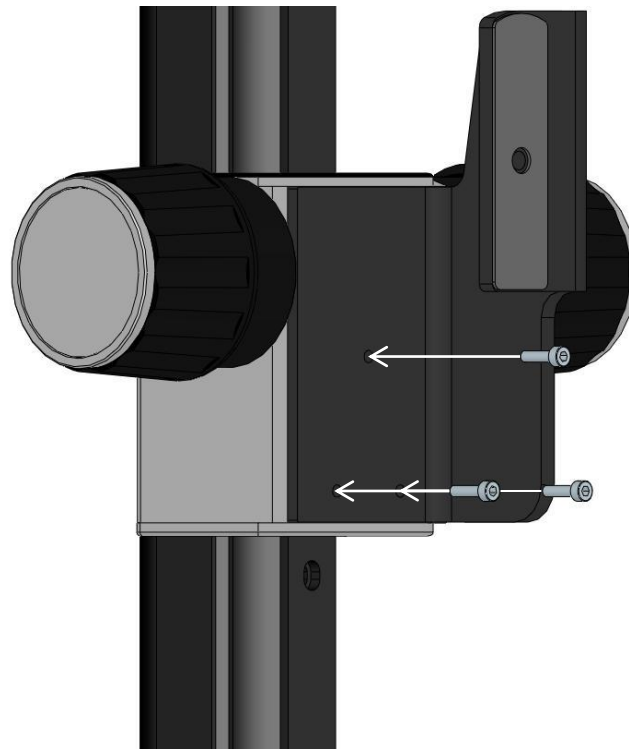

3. Slide in the Black and White Plate in the middle of the Base by pushing the spring on side.

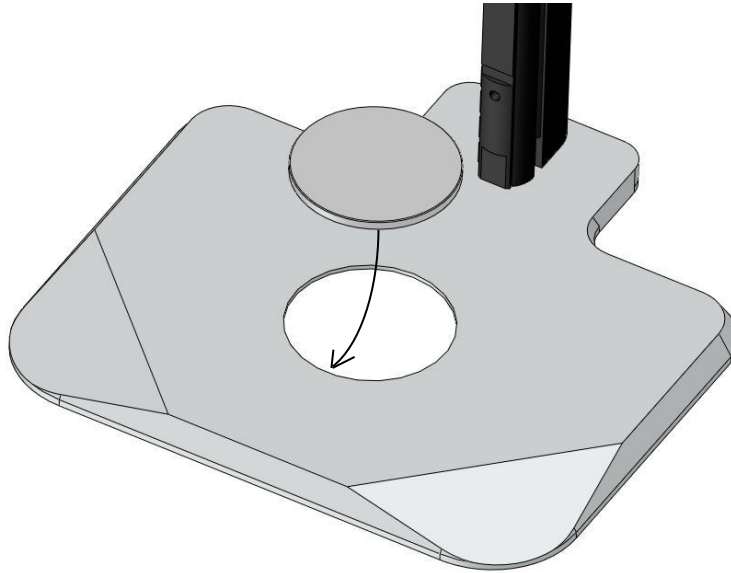

4. Cover the top with the Cover Cap.
5. Put the Focusing Stand on an even table after installation. Install the camera onto the adapter, and fasten the screw.

Notes

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

All rights reserved. The information contained herein is designed only for use with Optilia Instruments AB's Focusing Stand. Optilia Instruments AB is not responsible for any use of this information for other purposes. Optilia Instruments AB shall not be liable to the purchaser of this product or third parties for damages, losses, costs or expenses incurred by the purchaser or third parties as result of: accident, misuses, or abuse of this product or unauthorized modifications, repairs or alternations to this product, or failure to strictly comply with Optilia Instruments AB's operating and maintenance instructions. Optilia Instruments AB shall not be liable for any damages or problems arising from the use of any options or any consumable products other than those designated as original Optilia products or Optilia approved products by Optilia Instruments AB. Optilia Instruments AB shall not be held liable for any damages resulting from electromagnetic interference that occurs from the use of any interface cables or devises other than those provided by or designated as Optilia-approved Products by Optilia Instruments AB.

Djupdalsvägen 22
S-192 51, Sollentuna
Sweden
Phone:
Fax:
E-mail:

+46 (0)8 35 33 60
+46 (0)8 35 33 63
Info@optilia.eu

